

# IPAA

**NON STERILE**

**Isopropanol 70% v/v  
blended with Water For Injection  
(WFI) quality water**



**USE INSTRUCTIONS:**

Non-Sterile InSpec IPA is designed for spraying, wiping and mopping applications. Hold spray approximately 15cm to 20cm from area to be treated. 5L versions, pour into an appropriate container for mopping. Apply to surface to ensure complete coverage. For screw cap bottles use the entire contents after opening.

**TECHNICAL DATA:**

Appearance: Clear liquid free from particles. Specific Gravity (20°C): 0.872 to 0.883. Alcohol proportion: 70% +/- 2% v/v.

**STORAGE:**

Store upright in original closed containers away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is available in the Material Safety Data Sheet (MSDS).

**SAFE HANDLING:**

Always wear gloves and goggles or face protection. Always read the label and MSDS before use.

**MATERIAL COMPATIBILITY:**

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical File.

**BENEFITS:**

- Low residue
- Rapid action
- Bactericidal and Fungicidal
- 3 months validated 'in-use' shelf life
- Contents protected using 'bag-in-bottle' presentation with Premium range
- All formats manufactured with WFI quality water
- Each bottle packed with two outer bags
- Ideal surface and aseptic transfer disinfectants
- Manufactured to GMP

**TECHNICAL USER INFORMATION:**

InSpec IPA is a non-sterile ready to use disinfectant. This is a 70% v/v alcohol solution in WFI quality water.

Non-Sterile InSpec IPA is manufactured to GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment.

The solution is filtered through a 0.2 micron filter.

InSpec IPA non-sterile is supplied with Certificates of Conformity and Analysis.

All formats of IPA have a stabilised 24 month shelf life.

**Microbiological Minimum Efficacy Testing Contact**

EN 1276 Bacteria	5 minutes	EN 13697 Bacteria	5 minutes
EN 1650 Fungi	15 minutes	EN 13697 Fungi	15 minutes

**InSpec™ IPA Non-sterile Product**

Product	Pack Size	Case Size	Product Code
InSpec™ IPA 1L Non-Sterile	1L trigger spray	6 x 1L	NSIPAWFI-1L
InSpec™ IPA 5L Non-Sterile	5L screw cap	2 x 5L	NSIPAWFI-5L
InSpec™ IPA 500ml Non-Sterile	500ml trigger spray	8 x 500ml	NSIPAWFI-500ML

